Gebrauchsanleitung | Operating manual | Mode d'em-
ploi | Instrucciones de manejo | 操作手册

Ladeständer

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BRAND GMBH + CO KG
Otto-Schott-Str. 25
97877 Wertheim (Germany)

T +49 9342 808 0
F +49 9342 808 98000
info@brand.de
www.brand.de

Do you need more operating manuals and translations?
Please refer to <http://www.brand.de/om> or use the following
Quick Response Code:

The original operating manual is in German. Other languages
are translations of the original operating manual.

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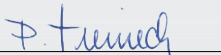
The manufacturer is solely responsible for issuing this declaration of conformity with regard to meeting the essential requirements and preparing the technical documentation.

| | |
|----------------------------------|--|
| Hersteller / Manufacturer | Adresse / Address |
| BRAND GMBH + CO KG | Otto-Schott-Str. 25 97877 Wertheim Germany |

Wertheim, 23. August 2022 / 23 August 2022

Patrick Ziemeck
 Managing Director Technology

i.A. Siegfried Ott
 Regulatory Affairs

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This document declares the accordance with the named harmonized regulations, but does not assure specific properties.

UK Declaration of Conformity

Product name: Charging stand for HandyStep® touch

Product variations: -

Cat.-No. 705220

| The product described above meets the basic requirements that are specified in the harmonization legislation listed below: | Applied harmonized standards: |
|---|---|
| The Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012 Radio Equipment Regulations 2017 (SI 2017 No. 1206, as amended by SI 2019 No. 696 Radio Equipment Regulations 2017 article 6.1a | EN IEC 63000:2018 EN 61010-1:2010/ A1:2019/AC:2019-04, EN 61010-1:2010/ A1:2019 |
| Radio Equipment Regulations 2017 article 6.1a Radio Equipment Regulations 2017 article 6.1b Radio Equipment Regulations 2017 article 6.1b Radio Equipment Regulations 2017 article 6.1b Radio Equipment Regulations 2017 article 6.2 | EN 62311:2008 EN 301 489-1 V2.1.1 EN 301 489-3 V2.1.1 EN 61326-1:2013 EN 303 417 V1.1.1 |
| The Approved Body: Element Materials Technology Warwick Ltd, Rothwell Road, Warwick, Warwickshire, CV34 5JX, UK AB number: 0891; performed a conformity assessment according Schedule 3, Module B and issued the RER Certificate: EMA21RER0035. Universal power supply not covered by the defined cert. | |
| Other applied standards: | Scope: |
| EN 62368-1:2014/AC:2015 (applied for power supply) 47 CFR §§ 18.305 / 18.307 KDB 680106 RSS-216 | LVD FCC (EMC USA) FCC (EMC USA) ISED (EMC Canada) |

The manufacturer is solely responsible for issuing this declaration of conformity with regard to meeting the essential requirements and preparing the technical documentation.

| Manufacturer | Address |
|--------------------|--|
| BRAND GMBH + CO KG | Otto-Schott-Str. 25 97877 Wertheim Germany |

Wertheim, 23 August 2022

Patrick Ziemeck
Managing Director Technology

i.A. Siegfried Ott
Regulatory Affairs

This document declares the accordance with the named harmonized regulations, but does not assure specific properties.

Declaration of conformity – China RoHS 2

BRAND GMBH + CO KG has made reasonable efforts to ensure that hazardous materials and substances may not be used in BRAND products.

In order to determine the concentration of hazardous substances in all homogeneous materials of the subassemblies, a “Product Conformity Assessment” (PCA) procedure was performed. As defined in GB/T 26572 the “Maximum Concentration Value” limits (MCV) apply to these restricted substances:

| | | | |
|-----------------|-------|---|------|
| + Lead (Pb): | 0.1% | + Hexavalent chromium (Cr(+VI)): | 0.1% |
| + Mercury (Hg): | 0.1% | + Polybrominated biphenyls (PBB): | 0.1% |
| + Cadmium (Cd): | 0.01% | + Polybrominated diphenyl ether (PBDE): | 0.1% |

Environmental Friendly Use Period (EFUP)

EFUP defines the period in years during which the hazardous substances contained in electrical and electronic products will not leak or mutate under normal operating conditions. During normal use by the user such electrical and electronic products will not result in serious environmental pollution, cause serious bodily injury or damage to the user's assets.

The Environmental Friendly Use Period for BRAND instruments is 40 years.

Material Content Declaration for BRAND Products

| 部件名称 Part name | 有毒有害物质或元素 Hazardous substances | | | | | |
|--|-----------------------------------|---------|---------|----------------|-------------|---------------|
| | 铅 Pb | 汞 Hg | 镉 Cd | 六价铬 Cr(+VI) | 多溴联苯 PBB | 多溴二苯醚 PBDE |
| 包装 / Packaging | 0 | 0 | 0 | 0 | 0 | 0 |
| 塑料外壳 / 组件 Plastic housing / parts | 0 | 0 | 0 | 0 | 0 | 0 |
| 电池 / Battery | 0 | 0 | 0 | 0 | 0 | 0 |
| 玻璃 / Glass | 0 | 0 | 0 | 0 | 0 | 0 |
| 电子电气组件 Electrical and electronic parts | X | 0 | 0 | 0 | 0 | 0 |
| 金属外壳 / 组件 Metal housing / parts | X | 0 | 0 | 0 | 0 | 0 |
| 电机 / Motor | X | 0 | 0 | 0 | 0 | 0 |
| 配件 / Accessories | X | 0 | 0 | 0 | 0 | 0 |
| 此表格是按照SJ/T 11364-2014中规定所制定的。 This table is created according to SJ/T 11364-2014. | | | | | | |

- O: 表示该有毒有害物质在该部件所有均质材料中的含量均在GB/T 26572规定的限量要求以下。
O: Indicates that the above mentioned hazardous substance contained in all homogeneous materials of the part is below the required limit as defined in GB/T 26572.
- X: 表示该有毒有害物质至少在该部件某一均质材料中的含量超出GB/T 26572规定的限量要求。
X: Indicates that the above mentioned hazardous substance contained in at least one of the homogeneous materials of this part is above the required limit as defined in GB/T 26572.

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Note: Batteries, glassware and accessories might not be content of the enclosed device and/or may have their own EFUP-marking and/or might be maintaining parts with changing EFUP-marking.

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Apart from the disclosures in the above table, the subassemblies are not intentionally manufactured or formulated with lead (Pb), mercury (Hg), cadmium (Cd), hexavalent chromium (Cr+VI), polybrominated biphenyls (PBB), and polybrominated diphenyl ethers (PBDE).

Products manufactured by BRAND may enter into further devices or can be used together with other appliances. With these third party products and appliances in particular, please note the EFUP labeled on these products. BRAND will not take responsibility for the EFUP of those products and appliances.

Place, date: Wertheim, 05|04|2022

Patrick Ziemeck
(Managing Director Technology)

i.A. Siegfried Ott
(Regulatory Affairs)

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Contents

| | | |
|----------|---------------------------------------|-----------|
| 1 | Scope of delivery | 16 |
| 2 | Safety | 16 |
| 3 | Function and Operation | 17 |
| 4 | Usage | 17 |
| 5 | Troubleshooting | 18 |
| 6 | Technical data | 18 |
| 6.1 | Charging Stand | 18 |
| 6.2 | Markings on Product and Battery | 18 |
| 7 | Order Information | 19 |
| 8 | Warranty | 20 |
| 9 | Disposal | 20 |

1 Scope of delivery

Charging stand (Use only with HandyStep® power supply/USB cable approved by BRAND).

2 Safety

Persons with medical implants must consult a physician before use in order to prevent the magnetic field emitted from influencing the implant.

Parts of the product become magnetic as soon as the power supply is established. Do not place magnetic data carriers (e.g. credit cards) near the charging stand.

Do not operate the charging stand outdoors and do not cover it.

USA

This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15B (HandyStep® touch and HandyStep® touch S) and part 18 (charging stand) of the FCC Rules.

These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications.

Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

Changes or modifications to this equipment not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

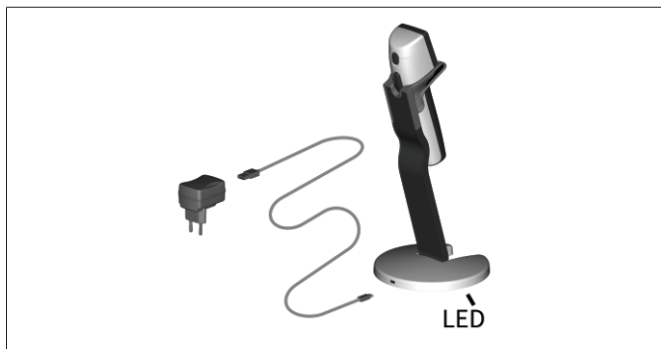
Canada

This device complies with the Industry Canada RSS-216 standard and operation is subject to the following two conditions:

1. This device may not cause interference.
2. This device must accept any interference received, including interference that may cause undesired operation of the device.

3 Function and Operation

The charging stand charges the HandyStep® inductively as soon as it has been placed in the charging stand.

4 Usage

Place HandyStep® in the charging stand as shown in the illustration and connect the power supply unit.

5 Troubleshooting

| Status LED | Device |
|------------------------|---|
| The LED's are lit. | HandyStep® is charging. |
| The LED's are not lit. | HandyStep® is not charging (battery charged or no power supply). |
| The LED's are flashing | An error has occurred. Charging is not possible. Reposition HandyStep® in charging stand. |

6 Technical data

6.1 Charging Stand

| | |
|---------------------|---|
| Input | DC 5.0 V  1.4 A, 7.0 W |
| Transmission | 3.5 W, 110 kHz ... 205 kHz |

6.2 Markings on Product and Battery

| Marking | Meaning |
|---|--|
|  | With this mark, we confirm that the product complies with the requirements set out in the EC Directives and has been subjected to the specified testing procedures. |
|  | Observe the information provided on the device, the accessories and in the operating instructions. |
|  (here: 40 years) | China RoHS (EFUP) EFUP defines the period in years during which the hazardous substances contained in electrical and electronic equipment do not leak or change under normal operating conditions. Under normal use by the user, such electrical and electronic products do not cause severe environmental pollution, serious personal injury or damage to the user's property. |
|  | DC voltage |

| Marking | Meaning |
|--|---|
|  | The electrical device must not be disposed of with the household waste. |
|  | UKCA: United Kingdom Conformity Assessed With this mark we confirm that the product complies with the requirements specified in the UK Designated Standards. |
| 2ATKA-HSTC705220 | Certification number USA FCC |
| IC: 25139-HSTC705220 HVIN: HW_HSTC01.00 | Certification number Canada ISED |

7 Order Information

| Designation | Figure | Ordering No. |
|--|---|--------------|
| Charging stand (worldwide except India) |  | 705220 |
| Support stand |  | 705230 |
| Universal AC Adapter HandyStep® touch and HandyStep® touch S incl. charging cable and country adapters |  | 705250 |

8 Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operating or unauthorized repairs of the device or for the consequences of normal wear and tear, especially of wearing parts such as pistons, seals, valves and the breakage of glass. The same applies for failure to follow the instructions of the operating manual. We are not liable for damage resulting from disassembly beyond that described in the operating manual or if non-original spare parts or components have been installed.

USA and Canada:

Find more warranty information on www.brandtech.com.

9 Disposal

Do not dispose of electrical appliances with household waste. Ensure that electrical equipment is disposed of properly.

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